

antibodies<sup>-online™</sup>

ABIN5311506

GST-Catcher

High-affinity anti-GST

Single-Domain Antibody (sdAb)

Protocol

---

For research use only

Not for use in clinical diagnostic procedures

Version Mar 2024

Catcher Product Line

GFP-Catcher - ABIN5311508

GFP-Catcher - ABIN7272855 Magnetic Beads

RFP-Catcher - ABIN5311510

RFP-Catcher - ABIN7529450 Magnetic Beads

BFP-Catcher - ABIN5311512

GST-Catcher - ABIN5311506

MBP-Catcher - ABIN7272855

mNeonGreen-Catcher - ABIN7529451

# Step-by-step Protocol

## I. Cell Collection & Lysis

1. For mammalian cells, harvest  $10^6$ - $10^8$  cells per sample.
2. Lyse cells according to established protocols in 0.2 to 1.5 mL volume.

Buffer recommendations:

2% Triton X-100, 1% Tween-20, 1% NP-40, 1% CHAPS, 1% Deoxycholate, 0.1% SDS


4 M NaCl, 2 M KCl, 1 M MgCl<sub>2</sub>, 100 mM EDTA

4 M urea

10 mM DTT, 10 mM 2-Mercaptoethanol


RNAse A, DNAse I, Benzonase, protease inhibitors

3. Centrifuge cell lysates in microcentrifuge tubes for 10 min at 14,000 x g
4. Transfer the supernatant to a fresh microcentrifuge tube for each sample and keep at 4 °C.

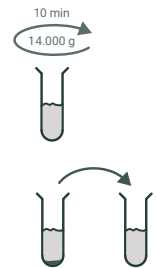


## II. Bead Preparation for GST Capture

5. Homogenize the GST-Catcher (agarose beads) slurry gently by shaking.
6. Transfer 20  $\mu$ L bead slurry to a 1.5 mL microcentrifuge tube for each sample.
7. Add 1 mL Lysis Buffer to equilibrate GST-Catcher (agarose beads).
8. Centrifuge GST-Catcher (agarose beads) for 1 min at 1000 x g and carefully remove the supernatant.
9. Repeat wash steps once for a total of two washes.

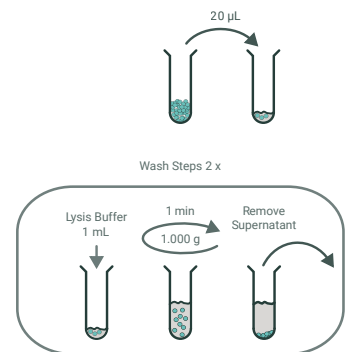

## III. Bead Incubation with Supernatant


10. Resuspend equilibrated GST-Catcher (agarose beads) gently with the cell lysate supernatant.
11. Rotate the microcentrifuge tubes for 1 h at 4 °C.
12. Centrifuge microcentrifuge tubes for 1 min at 1000 x g at 4 °C. Keep a small sample as "unbound" fraction. Carefully remove the supernatant.
13. Resuspend GFP-Catcher (agarose beads) in 1 mL Lysis Buffer.
14. Centrifuge GFP-Catcher (agarose beads) for 1 min at 1000 x g and carefully remove the supernatant.
15. Repeat wash steps twice for a total of three washes.

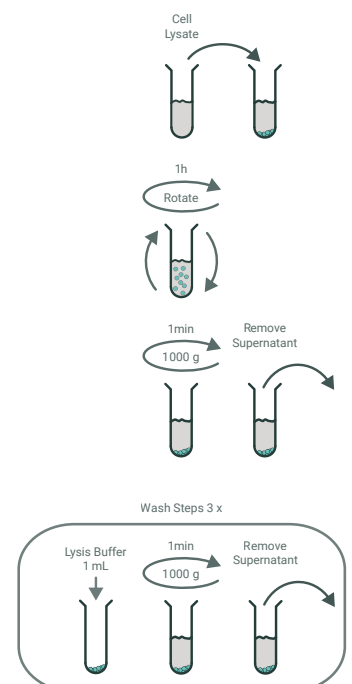

# Step-by-step Protocol

## IV. Bead Washing and Solution Changes

16. Resuspend GST-Catcher (agarose beads) gently in 1 mL TBS.
17. Centrifuge GST-Catcher (agarose beads) for 1 min at 1000 x g and carefully remove the supernatant.
18. Resuspend GST-Catcher (agarose beads) gently in 1 mL TBS.
19. Centrifuge GST-Catcher (agarose beads) for 1 min at 3000 x g and carefully remove the supernatant.

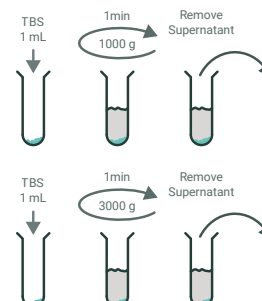

## V. Elution Preparation

20. Resuspend GST-Catcher (agarose beads) resin in 50  $\mu$ L 2X SDS sample buffer.
21. Heat sample (agarose beads) resin for 5 min to 95 °C.
22. Centrifuge microcentrifuge tubes for 1 min at 3000 x g and transfer the supernatant to fresh microcentrifuge tubes. Keep the pellet (agarose beads) as backup.



## For more information, please contact:

### antibodies-online Inc.

PO Box 5201  
Limerick, PA 19468  
USA

**Website:** [www.antibodies-online.com](http://www.antibodies-online.com)  
**Email:** [info@antibodies-online.com](mailto:info@antibodies-online.com)  
**Phone:** +1 877 302 8632  
**Fax:** +1 888 205 9894

### antibodies-online GmbH

Schloss-Rahe-Straße 15  
52072 Aachen  
Deutschland

**Website:** [www.antikoerper-online.de](http://www.antikoerper-online.de)  
**Email:** [info@antikoerper-online.de](mailto:info@antikoerper-online.de)  
**Phone:** +49 (0)241 95 163 153  
**Fax:** +49 (0)241 95 163 155